

Last Drop™ Low Retention Tips



SAMPLE RECOVERY
NOW BECOMES
EFFORTLESS


Features

- Manufactured from pure USP class VI certified virgin polypropylene free from natural rubber and inhibitory plasticizers
- 10 times less sample retention as compared to the standard pipette tips
- Full traceability
- Highly homogenous surface that prevents non-specific binding
- Diamond polished mould for the best-in-class hydrophobicity property
- Refill pack available in low retention universal pipette tips
- Available in all volume variants (Universal Pipette Tips and Universal Filter Pipette Tips), from 10 µl to 1250µl
- Colour coded inserts in racked tip box have details that include product code, product & volume description with lot number. This keeps your work organised by visual identification

Material & Method

- 200µl Last Drop™ pipette tips were compared with other low retention tips in combination with an electronic pipette, were used to aspirate & dispense 100µl of Malachite Green, Methylene Blue, Glycerol & Titron-X (liquid detergent) at a constant pipetting speed.
- Residual in the pipette were washed out by repeated rinsing with 200µl water into cuvette and compared photometrically. The average values established are shown in the table.

COMPARISON CHART

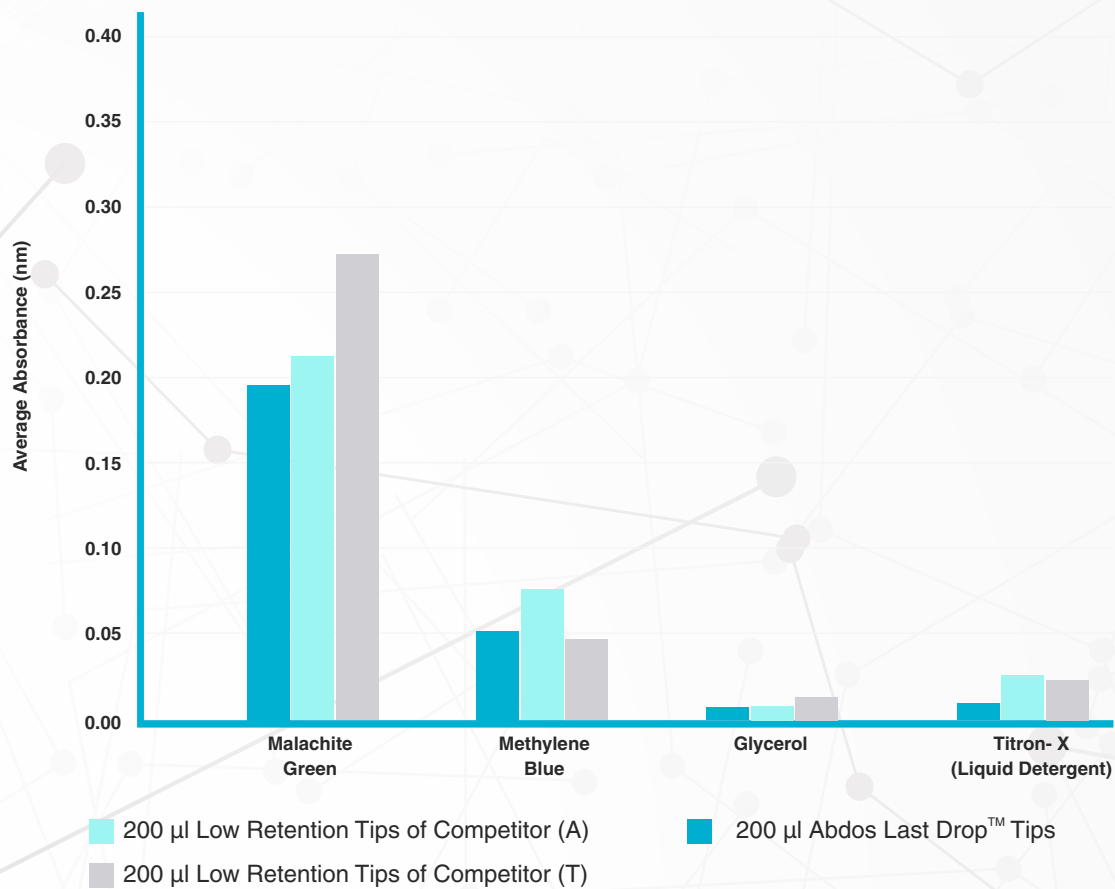

Result: Abdos Last Drop™ Pipette Tips clearly reduce sample loss. As seen above in the graph, it clearly indicates that Abdos Tips retained least amount of residue and saves on costly reagents.

Quality & Compliance



FREE



FREE



FREE



FREE



FREE




121°C



CERTIFIED

Bulk

- Easy to open and reseal
- Fully autoclavable
- Available in non-sterile

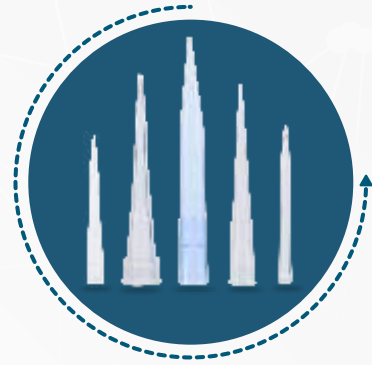

Individually Wrapped

- Fully automated blister packaging
- Gamma irradiated with SAL 10^{-6}
- Complete traceability with mention of product code, lot number and expiry date



Refill

- Saves precious time and storage space
- Refill inserts can be easily changed
- Available only in the non-filter category



Racked Tip Box

- Available sterile and non-sterile
- Colour coded rack inserts
- Design patented generation 2 tip box for stability
- Full traceability



Product Code	Product Description	No. per case
P10001	Last Drop™ Low Retention Pipette Tips, Bulk in Resealable Bags, 0.2 - 10 µl (Graduated Low Retention) Natural	1000
P10002	Last Drop™ Low Retention Pipette Tips, Bulk in Resealable Bags, 0.2 - 10 µl Extended (Graduated Low Retention) Natural	1000
P10003	Last Drop™ Low Retention Pipette Tips, Bulk in Resealable Bags, 2 - 200 µl (Bevelled Graduated Low Retention) Natural	1000
P10004	Last Drop™ Low Retention Pipette Tips, Bulk in Resealable Bags, 2 - 200 µl (Ultra Point Tip Orifice with Low Retention) Natural	1000
P10005	Last Drop™ Low Retention Pipette Tips, Bulk in Resealable Bags, 2 - 300 µl (Graduated Low Retention) Natural	1000
P10006	Last Drop™ Low Retention Pipette Tips, Bulk in Resealable Bags, 100 - 1000 µl (Low Retention) Natural	500
P10007	Last Drop™ Pipette Tips, Natural, 100 - 1000 XL (Extended)	500
P10008	Last Drop™ Pipette Tips, Natural, 100-1250, Last Drop Pipette Tips	500
P10010	Pure Right Last Drop™ Micro Tips, Individually Wrapped, Sterile, Natural, 0.2 - 10 µl (Graduated)	350
P10011	Pure Right Last Drop™ Micro Tips, Individually Wrapped, Sterile, Natural, 0.2 - 10 µl Extended (Graduated)	350
P10012	Pure Right Last Drop™ Micro Tips, Individually Wrapped, Sterile, Natural, 2- 200 µl (Bevelled Graduated)	350
P10013	Pure Right Last Drop™ Micro Tips, Individually Wrapped, Sterile, Natural, 2 - 300 µl (Graduated)	350
P10014	Pure Right Last Drop™ Micro Tips, Individually Wrapped, Sterile, Natural, 100 - 1000 µl	350
P10015	Pure Right Last Drop™ Micro Tips, Individually Wrapped, Sterile, Natural, 100 - 1000 µl XL (Graduated)	100
P10016	Pure Right Last Drop™ Micro Tips, Individually Wrapped, Sterile, Natural, 100 - 1250 µl (Graduated)	100
P10020	Last Drop™ Racked Micro, Tips 0.2 – 10 µl (Graduated Low Retention) Natural, 10 Racks x 96 Pcs	960

P10021	Last Drop™ Racked Micro, Tips 0.2 – 10 µl Extended (Graduated Low Retention) Natural, 10 Racks x 96 Pcs	960
P10022	Last Drop™ Racked Micro Tips, 2-200 µl (Bevelled Graduated Low Retention) Natural, 10 Racks x 96 Pcs	960
P10023	Last Drop™ Racked Micro Tips, 2-200 µl (Ultra Point Tip Orifice with Low Retention) Natural, 10 Racks x 96 Pcs	960
P10024	Last Drop™ Racked Micro Tips, 2-300 µl (Graduated Low Retention) Natural, 10 Racks x 96 Pcs	960
P10025	Last Drop™ Racked Micro Tips, 100-1000 µl (Low Retention) Natural, 10 Racks x 96 Pcs	960
P10026	Last Drop™ Racked Micro Tips, 100-1000 µl XL (Low Retention) Natural, 10 Racks x 96 Pcs	960
P10027	Last Drop™ Racked Micro Tips, 100-1250 µl (Low Retention) Natural, 10 Racks x 96 Pcs	960
P10040	Last Drop™ Racked Micro Tips 0.2 – 10 µl Sterile (Graduated Low Retention) Sterile Natural, 10 Racks x 96 Pcs	960
P10041	Last Drop™ Racked Micro, Tips 0.2 – 10 µl Extended (Graduated Low Retention) Sterile Natural, 10 Racks x 96 Pcs	960
P10042	Last Drop™ Racked Micro, Tips 2 - 200 µl (Bevelled Graduated Low Retention) Sterile Natural, 10 Racks x 96 Pcs	960
P10043	Last Drop™ Racked Micro Tips, 2-200 µl (Ultra Point Tip Orifice with Low Retention) Sterile Natural, 10 Racks x 96 Pcs	960
P10044	Last Drop™ Racked Micro Tips, 2-300 µl (Graduated Low Retention) Sterile Natural, 10 Racks x 96 Pcs	960
P10045	Last Drop™ Racked Micro Tips, 100 - 1000 µl (Low Retention) Sterile Natural, 10 Racks x 96 Pcs	960
P10046	Last Drop™ Racked Micro Tips, 100 - 1000 µl XL (Low Retention) Sterile Natural, 10 Racks x 96 Pcs	960
P10047	Last Drop™ Racked Micro Tips, 100 - 1250 µl (Low Retention) Sterile Natural, 10 Racks X 96 Pcs	960
P10060	Last Drop™ Refill Micro, Tips 0.2 - 10 µl (Graduated Low Retention) Natural, 12 Refill Racks x 96 Pcs	1152
P10061	Last Drop™ Refill Micro Tips 2-200 µl (Bevelled Graduated Low retention) Natural, 10 Refill Racks x 96 Pcs	960

P10062	Last Drop™ Refill Micro Tips 2-200 µl (Ultra Point Tip Orifice with Low Retention) Natural, 10 Refill Racks x 96 Pcs	960
P10063	Last Drop™ Refill Micro Tips 2-300 µl (Graduated Low Retention) Natural, 10 Refill Racks x 96 Pcs	960
P10064	Last Drop™ Micro Tips – Refill Pack, PP, 100 - 1000 µl (XL), 6 Refill Racks x 96 Pcs	576
P10065	Last Drop™ Micro Tips – Refill Pack, PP, 100 - 1250 µl Natural, 6 Refill Racks x 96 Pcs	576
P11050	Last Drop™, Low Retention, Filter Micro Tips, PP, 0.2 - 10 µl (Natural) (Graduated), Bulk	1000
P11051	Last Drop™, Low Retention, Filter Micro Tips, PP, 0.2 - 10 µl Extended (Natural) (Graduated), Bulk	1000
P11052	Last Drop™, Low Retention, Filter Micro Tips, PP, 2-20 µl (Natural) (Bevelled Graduated), Bulk	1000
P11053	Last Drop™, Low Retention, Filter Micro Tips, PP, 2-100 µl (Natural) (Bevelled Graduated), Bulk	1000
P11054	Last Drop™, Low Retention, Filter Micro Tips, PP, 2-200 µl (Natural) (Bevelled Graduated), Bulk	1000
P11055	Last Drop™, Low Retention, Filter Micro Tips, PP, 2-300 µl (Natural) (Graduated), Bulk	1000
P11056	Last Drop™, Low Retention, Filter Micro Tips, PP, 100-1000 µl (Natural), Bulk	500
P11057	Last Drop™, Last Drop Filter Pipette Tips, PP, 100 - 1000 µl, XL (Extended)	500
P11058	Last Drop™, Last Drop Filter Pipette Tips, PP, 100 - 1250 µl	500
P11040	Pure Right Last Drop™ Low Retention Filter Tips; Individually Wrapped, 10 ul	350
P11041	Pure Right Last Drop™ Low Retention Filter Tips; Individually Wrapped, 10 ul, Extended	350
P11042	Pure Right Last Drop™ Low Retention Filter Tips; Individually Wrapped, 20 ul	350
P11043	Pure Right Last Drop™ Low Retention Filter Tips; Individually Wrapped, 100 ul	350

P11044	Pure Right Last Drop™ Low Retention Filter Tips; Individually Wrapped, 200 ul	350
P11045	Pure Right Last Drop™ Low Retention Filter Tips; Individually Wrapped, 300 ul	350
P11046	Pure Right Last Drop™ Low Retention Filter Tips; Individually Wrapped, 1000 ul	350
P11070	Pure Right Last Drop™ Low Retention Filter Tips; Individually Wrapped, 1000 ul, XL	100
P11071	Pure Right Last Drop™ Low Retention Filter Tips; Individually Wrapped, 1250 ul	100
P11060	Last Drop™, Low Retention, Racked Filter Micro Tips, 0.2-10 µl (Natural) (10 Racks X 96 Pcs)	960
P11062	Last Drop, Low Retention, Racked Filter Micro Tips, 0.2-10 µl Extended (Natural) (10 Racks X 96 Pcs)	960
P11064	Last Drop™, Low Retention, Racked Filter Micro Tips, 2-20 µl (Natural) (Bevelled Gradated) (10 Racks X 96 Pcs)	960
P11066	Last Drop™, Low Retention, Racked Filter Micro Tips, 2-50 µl (Natural) (Bevelled Gradated) (10 Racks X 96 Pcs)	960
P11068	Last Drop™, Low Retention, Racked Filter Micro Tips, 2-100 µl (Natural) (Bevelled Gradated) (10 Racks X 96 Pcs)	960
P11072	Last Drop, Low Retention, Racked Filter Micro Tips, 2-200 µl (Natural) (Bevelled Gradated) (10 Racks X 96 Pcs)	960
P11074	Last Drop™, Low Retention, Racked Filter Micro Tips, 2-300 µl (Natural) (Gradated) (10 Racks X 96 Pcs)	960
P11076	Last Drop™, Low Retention, Racked Filter Micro Tips, 100-1000 µl (Natural) (10 Racks X 96 Pcs)	960
P11078	Last Drop™, Low Retention, Racked Filter Micro Tips, 100-1000 XL, Extended, (Natural) (10 Racks X 96 Pcs)	960
P11080	Last Drop™, Low Retention, Racked Filter Micro Tips, 100-1250 µl (Natural) (10 Racks X 96 Pcs)	960
P11061	Last Drop™, Low Retention, Racked Filter Micro Tips, 0.2-10 µl (Natural), Sterile (Gama Radiation) (10 Racks X 96 Pcs)	960
P11063	Last Drop™, Low Retention, Racked Filter Micro Tips, 0.2-10 µl Extended (Natural), Sterile (Gama Radiation) (10 Racks X 96 Pcs)	960

P11065	Last Drop™, Low Retention, Racked Filter Micro Tips, 2-20 µl (Natural), Sterile (Bevelled Gradated) (Gama Radiation) (10 Racks X 96 Pcs)	960
P11067	Last Drop™, Low Retention, Racked Filter Micro Tips, 2-50 µl (Natural), Sterile (Bevelled Gradated) (Gama Radiation) (10 Racks X 96 Pcs)	960
P11069	Last Drop™, Low Retention, Racked Filter Micro Tips, 2-100 µl (Natural), Sterile (Bevelled Gradated) (Gama Radiation) (10 Racks X 96 Pcs)	960
P11073	Last Drop™, Low Retention, Racked Filter Micro Tips, 2-200 µl (Natural), Sterile (Bevelled Gradated) (Gama Radiation) (10 Racks X 96 Pcs)	960
P11075	Last Drop™, Low Retention, Racked Filter Micro Tips, 2-300 µl (Natural), Sterile (Gradated) (Gama Radiation) (10 Racks X 96 Pcs)	960
P11077	Last Drop™, Low Retention, Racked Filter Micro Tips, 100 -1000 µl (Natural), Sterile (Gama Radiation), (10 Racks X 96 Pcs)	960
P11079	Last Drop™, Low Retention, Racked Filter Micro Tips, 100 - 1000 XL (Extended), (Natural), Sterile (Gama Radiation) (10 Racks X 96 Pcs)	960
P11081	Last Drop™, Low Retention, Racked Filter Micro Tips, 100 -1250 µl (Natural), Sterile (Gama Radiation) (10 Racks X 96 Pcs)	960

What Abdos Life Sciences stands for?



Precision



Reliability



Responsiveness



Innovation



Sustainability



Relationships

Get in touch with us:

📞 011-41081897

✉️ labtech@abdosindia.com

🌐 www.abdoslifesciences.com