

Technical Data Sheet

Doc No– Ab/TDS/Eqpt -233

Rev No : 05 Date 02/2022-23

TECHNICAL DATA SHEET

Abdos Labtech Pvt. Ltd.

SWIRLTOP – 7inch LED Digital Magnetic Stirrer & Hot Plate

Valid for Catalog No. E11650, E11652, E11651, E11653

1. Description/Specification

1.1 Description: SWIRLTOP – 7inch LED Digital Magnetic Stirrer & Hot Plate

1.1.1 LED display for precise monitoring of speed and temperature for Magnetic stirrer & Hot plate.

1.1.2 Built in PID Controller ensures the safe heating of the medium with overheating and protection. Max heating plate temperature 550°C.

1.1.3 Glass ceramic work plate provides excellent chemical resistant performance and efficient heat transfer.

1.1.4 External temperature control is possible by connecting the temperature sensor (PT 1000) with an accuracy at $\pm 0.2^{\circ}\text{C}$.

1.1.5 Digital speed control. Maximum Speed at 1500rpm. Outstanding brushless DC motor performance enables more stirring power.

1.1.6 Two rotating knobs enable the easy and direct change of speed and temperature.

1.1.7 HOT WARNING will flash when the work plate temperature is above 50°C or even the hot plate is turned off.

1.1.8 Remote Function provides PC Control and data transmission via RS232 using 3rd Party Software.

1.2 Design: SWIRLTOP – 7inch LED Digital Magnetic Stirrer & Hot Plate

1.3 Performance Specifications:

Specifications	E11650 & E11652	E11651 & E11653
Voltage	220-240V	220-240V
Frequency	50-60Hz	50-60Hz
Over Heating Protection	580°C	580°C
Speed Range	100-1500 rpm	100-1500 rpm
Overload Protection	LED Lights, motor automatically stops	LED Lights, motor automatically stops
Motor Protection	LED Lights, motor automatically stops	LED Lights, motor automatically stops
Temp. Display	LED	LED
Speed Display	LED	NA
Max. Stirrer Quantity	20L	NA



Abdos Labtech Private Limited
(A Division of Abdos Group)
(An ISO 9001,13485,14001,18385,GMP & CE certified company)
Shiv Ganga Industrial Estate
Plot No-1, Lakheswari, Bhagwanpur,
Roorkee-247667, (UK) INDIA

Registered office
13A Government
Place East, 2nd Floor Kolkata
West Bengal – 700069
CIN: U85100WB2008PTC123903
Corporate Office
319, 3rd Floor, DLF Tower A, DDA Dist. Centre
Jasola, New Delhi - 110025

Ph. No: +91- (0) 1332235277
Email: labtech@abdosindia.com
Web: www.abdoslivesciences.com
GST: 05AAGCA8605A1Z7

Technical Data Sheet

Max. Magnetic bar length	80mm	NA
Work Plate Material	Glass Ceramic	Glass Ceramic
Dimension	83x220x186 mm	83x220x186 mm
Permissible ambient Temperature	5 - 40°C	5 - 40°C
Permissible ambient	80%	80%
Protection class according	IP21	IP21

Code	Product	Workplace size (cm)
E11650	Magnetic Stirrer & Hot Plate with Support rod & external temp. probe	18.4x18.4
E11651	Hotplate with support rod & external temp. probe	18.4x18.4
E11652	Magnetic Stirrer & Hot Plate	18.4x18.4
E11653	Hot Plate	18.4x18.4

Spare Accessories:

Code	Product
E11654	SP-Support rod 7inch
E11233	External temp. probe

1.4 Material: Body – ABS Material

1.5 Color: Grey

1.6. Basic features:

1.6.1 Two rotating knobs enable the easy and direct change of speed & Temp.

1.6.2 Digital Speed Controlling.

1.6.3 Max. Heating plate temp. 550 °C

1.6.4 Remote function provides PC control and data transmission.

1.6.5 **Material:** Made up of highly pure Virgin ABS (Acrylonitrile Butadiene Styrene), Glass Ceramic.

1.6.6 CE certified

1.6.7 **Chemical resistance:** see website (www.abdoslifesciences.com)



Abdos Labtech Private Limited
(A Division of Abdos Group)
(An ISO 9001,13485,14001,18385,GMP & CE certified company)
Shiv Ganga Industrial Estate
Plot No-1, Lakheswari, Bhagwanpur,
Roorkee-247667, (UK) INDIA

Registered office
13A Government
Place East, 2nd Floor Kolkata
West Bengal – 700069
CIN: U85100WB2008PTC123903
Corporate Office
319, 3rd Floor, DLF Tower A, DDA Dist. Centre
Jasola, New Delhi - 110025

Ph. No: +91- (0) 1332235277
Email:labtech@abdosindia.com
Web: www.abdoslifesciences.com
GST: 05AAGCA8605A1Z7

Technical Data Sheet

Warranty: 12 months from the date of Installation or 15 months from the date of Invoicing

Instruction for Use

Storage temp. Max 30°C

Before using the product check crackness, pouch punctured etc.

Check all the information printed on pouches or boxes.

WARNING

DO NOT store strong oxidizing agents.

DO NOT place any plastic labware in a flame.

DO NOT mix any chemicals that may result in a thermic reaction, which can cause product failure

Consult local fire codes prior to storage of flammable liquids.



Abdos Labtech Pvt. Ltd.

Plot No. 1 Lakeshwari Bhagwanpur Roorkee 247667 (UK) India
Email : labtech@abdosindia.com, www.abdoslifesciences.com
Ph: 4(91) 11 41081897



Quality Assurance Incharge
Abdos Labtech Pvt. Ltd.



Abdos Labtech Private Limited
(A Division of Abdos Group)
(An ISO 9001,13485,14001,18385,GMP & CE certified company)
Shiv Ganga Industrial Estate
Plot No-1, Lakshewari, Bhagwanpur,
Roorkee-247667, (UK) INDIA

Registered office
13A Government
Place East, 2nd Floor Kolkata
West Bengal – 700069
CIN: U85100WB2008PTC123903

Corporate Office
319, 3rd Floor, DLF Tower A, DDA Dist. Centre
Jasola, New Delhi - 110025

Ph. No: +91- (0) 1332235277
Email: labtech@abdosindia.com
Web: www.abdoslifesciences.com
GST: 05AAGCA8605A1Z7

Technical Data Sheet



Abdos Labtech Private Limited
(A Division of Abdos Group)
(An ISO 9001,13485,14001,18385,GMP & CE certified company)
Shiv Ganga Industrial Estate
Plot No-1, Lakheswari, Bhagwanpur,
Roorkee-247667, (UK) INDIA

Registered office
13A Government
Place East, 2nd Floor Kolkata
West Bengal – 700069
CIN: U85100WB2008PTC123903

Corporate Office
319, 3rd Floor, DLF Tower A, DDA Dist. Centre
Jasola, New Delhi - 110025

Ph. No: +91- (0) 1332235277
Email:labtech@abdosindia.com
Web: www.abdoslivesciences.com
GST: 05AAGCA8605A1Z7